




**CAUTION:** This temperature sensor is sensitive to electrostatic discharge (ESD). Use ESD precautionary procedures when handling, or making mechanical or electrical connections to this device in order to avoid performance degradation or loss of functionality.

There are three aspects of using a temperature sensor which are critical to its optimum performance. The first involves the proper mounting of the sensor package, the second relates the proper joining of sensor lead wires and connecting wires. The final concern is the thermal anchoring of the lead wires. Although the sequence in which these areas should be addressed is not set in stone, all elements covered under each aspect should be adhered to for maximum operating capabilities of the sensor.

### SENSOR MOUNTING

The DT-470/471/670/671 LR package combines a standard SD sensor mounted on a gold plated copper adapter and is designed for mounting in a 3.2 mm (0.125 in) hole.

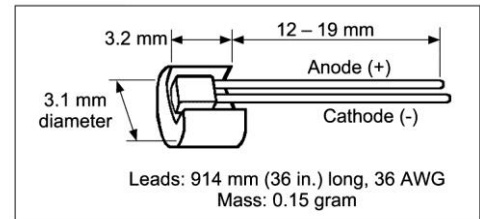

1. A hole should be drilled 3.2 mm (0.125 in) diameter by 3.2 mm (0.125 in) deep minimum for the LR adapter.
2. Surface area should be cleaned with a solvent such as Acetone followed by an Isopropyl Alcohol rinse. Allow time for the solvents to evaporate before sensor positioning.
3. A small amount of Apiezon<sup>®</sup> N grease should be applied around the mounting surface and the sensor to enhance thermal contact.

**Note:** It is important that no electrically conductive particles lodge between the adapter and sensor. There is a thin braze joint around the sides of the SD sensor and is electrically connected to the sensing element. Contact to the sides with any electrically conductive material will cause a short.


4. Position the sensor package so the LR adapter is fully submersed in the mounting hole.

**Note:** The LR package is designed for use up to 420 K (147 °C) with the standard curve, but the standard calibration only goes to 325 K (52 °C). Consult Lake Shore for calibrations above 325 K (52 °C).

**Note:** Be careful that the sensor leads do not touch the mounting surface. This will cause a short.

### LEAD ATTACHMENT

It is recommended that a 4-lead measurement scheme is used with this sensor. Attaching four connecting wires to the sensor leads is recommended. See figure to determine sensor lead polarity.

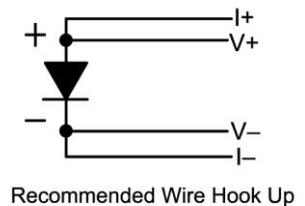

1. Prepare the sensor leads and connecting lead wires with a RMA (rosin mildly active) soldering flux, tin them with a minimal amount of 60% Sn 40% Pb solder. Use a low wattage soldering iron that will not exceed 200 °C. Clean off residual flux with rosin residue remover. The sensing element inside the package should be protected from excessive heat by putting a heat sink clip over the package.
2. Strip connecting wires insulation by delicately scraping with a razor blade, fine sand paper or steel wool. Phosphor-bronze or manganin wire, in sizes 32 or 36 AWG, is commonly used as the connecting lead wire. These wires have low thermal conductivity which help minimize the heat flow through the leads. Typical wire insulation is Polyvinyl Formal (Formvar<sup>™</sup>) or Polyimide (ML). Formvar<sup>™</sup> insulation has better mechanical properties such as abrasion resistance and flexibility. Polyimide insulation has better resistance to chemical solvents and burnout.
3. Prepare the connecting wire ends with a RMA (rosin mildly active) soldering flux, tin them with a minimal amount of 60% Sn 40% Pb solder. Use a low wattage soldering iron which will not exceed 200 °C.
4. Clean off residual flux with rosin residue remover. The sensor lead can be prepared in an identical manner.
5. Attach one sensor lead with two of the connector wires and apply the soldering iron above the joint area until the solders melt, then remove the iron immediately. Repeat for the other set of connector wires and the other sensor lead.
6. Avoid putting stress on the device leads and leave enough slack to allow for the thermal contractions that occur during cooling which could fracture a solder joint or lead. This can be achieved with heat shrink tubing.

## HEAT SINKING/THERMAL ANCHORING

1. Since the area being measured is read through the base of the sensor, heat flow through the connecting leads can create an offset between the sensor chip and the true sample temperature. Thermal anchoring of the connecting wires is necessary to assure that the sensor and the leads are at the same temperature as the sample.
2. Connecting wires should be thermally anchored at several temperatures between room temperature and cryogenic temperatures to guarantee that heat is not being conducted through the leads to the sensing element.
3. If the connecting leads have a thin insulation such as Formvar™ or Polyimide, a simple thermal anchor can be made by winding the wires around a copper post, bobbin or other thermal mass. A minimum of five wraps around the thermal mass should provide enough of an anchor. However, if space permits, additional wraps are recommended for good measure. To maintain good electrical isolation over many thermal cycles, it is good practice to first varnish a single layer of cigarette paper to the anchored area then wrap the wire around the paper and bond in place with a thin layer of IMI 7031 Varnish. Formvar™ wiring insulation has a tendency to craze with the application of IMI varnish. Once IMI varnish is applied, the wires cannot be disturbed until all solvents have evaporated and the varnish has fully cured (typically 12 to 24 h).
4. A final thermal anchor at the sample itself is a good practice to ensure thermal equilibrium between the sample and temperature sensor.

**CRYOGENIC ACCESSORIES** – Recommended for proper installation and use of DT-470/471/670/671 LR Sensors.

**Stycast® Epoxy 2850FT** (P/N 9003-020, 9003-021): Permanent attachment, excellent low temperature properties, poor electrical conductor, low cure shrinkage.

**Apiezon® N Grease** (P/N 9004-020): Low viscosity, easy to use, solidifies at cryogenic temperatures, excellent lubricant.

**IMI 7031 Varnish** (P/N 9009-002): Nonpermanent attachment, excellent thermal conductor, easy to apply and remove.

**Indium Solder** (P/N 9007-002-05): 99.99% pure, excellent electroplating material, foil form.

**90%Pb 10%Sn Solder** (P/N 9008-001): Greater lead content, for higher temperature applications no greater than 200°C.

**RMA Soldering Flux** (P/N 9008): Variety of types, refer to Lake Shore Product Catalog for details.

**Phosphor-bronze Wire** (P/N 9001-00X): Available in single, duo, and quad strands, no magnetic attraction, low thermal conduction.

**Manganin Wire** (P/N 9001-00X): Low thermal conductivity, high resistivity, no magnetic attraction.

**Heat Sink Bobbin** (P/N 9007-900 Large, 9007-901 Small): Gold plated oxygen-free high-conductivity (OFHC) copper bobbins.

**Instruments:** Lake Shore sells a complete line of instrumentation used with the DT-470/471/670/671 sensors, such as: current sources, cryopump monitors, temperature controllers, monitors and thermometers, temperature scanners and transmitters.

For complete product description and detailed specifications on the above accessories and instruments, consult the Lake Shore Temperature Measurement and Control Catalog, call at (614) 891-2243, e-mail at [sales@lakeshore.com](mailto:sales@lakeshore.com), or visit our website at [www.lakeshore.com](http://www.lakeshore.com).